


440-Z

440G X 0.1G CAPACITY

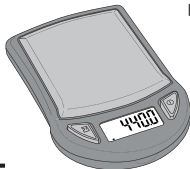

ENGLISH



USER MANUAL

Thank you for purchasing the My Weigh® 440-Z digital pocket scale. Please read all operating instructions carefully before use. This electronic scale is a precision instrument. With normal care and proper treatment, it will provide years of reliable service. For more information please visit www.myweigh.com.

Never load the scale with more than the maximum capacity. Although the 440-Z is designed to be extremely durable with extra overload protection built into the case, overloading will permanently damage it! Avoid any exposure to extreme heat or cold, your scale works better when operated at normal room temperature. Keep your scale in a clean environment. Dust, dirt, moisture, vibration, air currents and/or a close proximity to other electronic equipment can all cause an adverse effect on the reliability and accuracy of your scale. Handle with care. Gently apply all items to be weighed onto tray top. Avoid shaking, dropping or otherwise shocking the scale. Scales are delicate instruments and unlike cellular phones, scales have delicate sensors that determine how much an item weighs. If you drop or shock your scale, these sensors "feel" the shock and are sometimes destroyed. This happens with all digital scales. We design our scales to be as resistant to shock or drops as possible, however there is no way for us to protect 100% against load cell or sensor damage.

Failure to follow these instructions will void your warranty.

Always allow the scale to acclimate to normal room temperature for at least one hour before use. Give your scale sufficient warm up time. Usually 30-60 seconds before calibration to give the internal components a chance to stabilize.

BATTERIES

Low Batteries & bad battery connections are the #1 cause of scale malfunction and inaccuracy! We test **all** of our scale returns from consumers and 60% of them are battery related problems. This sounds **silly** but it's true! A scale **will** perform **poorly** when it has **low** batteries. Use **good quality** batteries & **replace** them often (Remove the batteries if you **plan** to store the scale for **longer** then 14 days). We **include** good quality batteries with **all** of our scales but they can run **low** in storage. If your scale simply won't turn on while on battery power, it is often caused by **loose** battery connections.



Battery installation

- Press and **lift** open the battery cover **located** at the bottom of the unit.
- Insert** the batteries and make sure the **polarity** is correct.
- Close** the battery cover until it **clicks** shut.



Note : If the battery symbol appears in the **display**, it means **low** battery power. It is time to **replace** the batteries.

OPERATION INSTRUCTIONS

Weighing Procedures

- Press [] to turn on or off the scale.
- Press [] to select a weighing unit g, oz, ozt, dwt.
- Start weighing.

Tare




Tare can be used for **eliminating** the weight **value** of an empty container. **Place** an empty container on the scale and press []. After 3 seconds **place** the items to be weighed in the container. NOTE: When **all** weight is removed from the weighing tray, the **tared** value of a container **will** be **displayed** as a negative number. Press [] again to return the scale to zero.

CALIBRATION

When to **calibrate** - calibration is recommended for best results.

Calibration may be required when the scale is first set up for use, or if the scale is moved to a different **altitude** or new **location**. This is necessary because the weight of a mass in one **location** is not necessarily the same in another **location**. Also, with time and use, mechanical **deviations** can occur.

To **Calibrate** the 440-Z: you must have an accurate 200g weight in order to **calibrate**.

1. Turn the scale **OFF**, slide the switch on the bottom of the scale from "Norm" to "CAL". Place the scale on a **Flat**, very **stable** surface (Be certain the scale is **OFF**).
2. Press the [] key the display will show the a/d value .
3. Press the [] key, the display will read "zero" and then the correct **calibration** weight
4. Gently place the 200g weight on the tray and wait for 3 seconds.
5. Press the [] key. The display will show "CAL", "PASS", and then will return to normal weighing and show 200.0
6. Calibration is now complete. Remove the weight and slide the switch on the bottom of the scale from "CAL" to "Norm".

FEATURES


Power Up Segment Test

When first turning the unit on, the scale will run a quick diagnostic and it will display a countdown . This display will remain for a few seconds and then reset to zero.

Overload indicator

When the display shows "EEEE", this indicates an **overload**. Remove excessive load immediately. Remember: you can permanently damage the scale and void your warranty by **overloading** it!

Negative Value

Any tared value will be displayed as a negative number once all weight is removed, press [] to re-zero the scale.

Auto off

An auto off feature is provided to conserve battery power. The unit will automatically turn off after 2 minutes of inactivity.

Expansion Tray/Cover

The protective cover of the 440-Z is designed to also be used as a platform expansion tray.

Capacity	440g x 0.1g	Units	g, oz, ozt, dwt
Auto-OFF	2 min.		
Scale dimension	63mm x 94mm x 20mm		
Tray dimension	53mm x 63mm		
Net/gross weight	54g		
Operating temperature	Optimum 10-30°C (50-86°F)		
Power Source	2 x AAA alkaline batteries		

